

TÍTULO: FLASH FROZEN PRESERVATION OF LUNG TISSUE IN OCT		
ARCHIVADO: PNT_8.2.006_v1.3_EN_LUNG_TISSUE_FLASH_FROZEN_OCT		Nº DE CÓDIGO: 8.2.006
VERSIÓN: 1.3	FECHA DE EDICIÓN: 26/07/2017 11:07:00 O7/P7	PÁG. 1 DE 4
REVISADO POR: Laura Vidaña, Estefanía Piñero, Victor Miguel Galvez, Daniel Pons FECHA: 24/07/2017		FIRMA:
APROBADO POR: Cristina Villena FECHA: 26/07/2017		FIRMA:

1. OBJECTIVE

Tissue samples are collected from surgical surplus from patients that have signed the informed consent form and agree with the participation in the Pulmonary Biobank Consortium.

Fresh frozen lung tissue collections are a valuable resource for research. Tissues suitably preserved are appropriate for proteomic and genomic studies. OCT (*Optimal Cutting Temperature*) is a specific compound used to preserve ultra-cellular structure (histology) and probably nucleic acids structure. This document aims to normalize and standardize lung tissue freezing in OCT protocol for Pulmonary Biobank Consortium.

2. SCOPE

This protocol describes how to quickly freeze lung tissue in OCT. This procedure does not detail biohazard and/or chemicals processes for occupational safety and health. It is recommended that the staff follows the established Health and Safety rules of each center.

3. RELATED DOCUMENTS

PNT_8.2.001_Recogida y transporte Tejido Pulmonar.doc
PNT_8.1.001_Etiquetado y Trazabilidad.doc

4. ROLES AND RESPONSIBILITIES

The fulfilment of this standard protocol rests with all the members of the Pulmonary Biobank Consortium who are responsible for the flash freezing of lung tissue in OCT.

5. MATERIALS AND EQUIPMENT

TÍTULO: FLASH FROZEN PRESERVATION OF LUNG TISSUE IN OCT**ARCHIVADO:**

PNT_8.2.006_v1.3_EN_LUNG_TISSUE_FLASH_FROZEN_OCT

Nº DE CÓDIGO: 8.2.006

Recommended materials and equipment are listed below. Depending on the place where the task or procedure is performed, these materials may be replaced by alternative or equivalent products.

Materials and equipment	Materials and equipment (Specific to Center)
Containers for liquid nitrogen transportation	
Liquid nitrogen	
2-methylbutane (isopentane)	
Isopentane container (bath in liquid nitrogen or dry ice)	Histobath or equivalent
Suitable containers for handling and cutting (Petri dishes)	
Pens and Markers	
Sterile forceps	
PBS or saline solution	
Vinyl cryomolds (such as Tissue-Tek Ref.#4557 25mm x 20 mm x 5 mm)	
Freezing medium OCT (Tissue-Tek O.C.T. Ref.#4583)	
Scissors to cut tissue	
Gloves to protect staff from tissue handling	
Adequate and suitable labels for cryomolds	
Splash protection mask	
Lab coat for protection against spills and splashes	
Protection mask suitable for liquid nitrogen handling	
Gloves suitable for liquid nitrogen handling	

6. GLOSSARY

Cryopreservation: Biological conservation process that consists on the preservation at very low temperatures for a long period of time.

OCT: "*Optimal Cutting Temperature*" refers to a compound used to preserve tissues against freezing. OCT preserves ultra-structures and protects the tissue from dehydration and degradation. It acts as an isolator on temperature variations and minimizes crystal formation. It is especially useful for preservation of frozen fresh tissue that will be sectioned.

7. PROCEDURE

The objective of this procedure is to guarantee safety, timely and efficient tissue sample preservation from patients that have signed the informed consent form and agree with the participation in the Pulmonary Biobank Consortium. Thereby this protocol tries to avoid contamination risk and molecular integrity loss. The sample preservation process is essential for

high integrity and quality products obtainment, which facilitate the use of genomic and proteomic techniques.

7.1 TISSUE TRANSPORTATION

1. Tissue mobilization must be quick; no more than 30 minutes between resection and the beginning of tissue conservation. To ensure it, manual or digital registration of time should be performed.
2. If time between resection and the beginning of tissue conservation is over 30 minutes, it should be indicated as “incidence” in the sample information field or in the “observations” field. If time exceeds 40 minutes, the tissue collection according to this protocol should not be done.
3. The operating room staff will notify the tissue availability to the person responsible of tissue obtainment who will organize the tissue transport from the operating room to the assigned lab. Transportation will be organized optimally to preserve cellular and molecular integrity.
4. The pathologist or responsible delegated person will determine the tissue fragment surplus for biobank based on clinical needs.
5. Tissue sample must be introduced into a plastic container, which in turn is inside a container filled with ice.
6. If possible, prepare the tissue collection kits previously.

7.2 FLASH FREEZING OF LUNG TISSUE IN OCT

1. Preservation is carried out by the laboratory technician or the person designated by the head of each Pulmonary Biobank Consortium node.
2. Treat the tissue as potentially infectious and do not place it directly on ice. Place tissue in PBS or saline solution pre-cooled at 4°C in a 50ml sterile container.
3. Wash 3 times with pre-cooled PBS or saline solution, shaking for 3-4 minutes or static for 10 minutes.
4. Cut the sample in approximately 0,3 x 0,3 x 0,3 cm fragments with the aid of sterile scissors and forceps.
5. Lightly dry the sample on filter paper.
6. Make sure that sectioned tissue does not dry out or become contaminated by contact with other tissues or samples. Use always clean and sterile forceps and scissors. Avoid cross contamination between samples or between tumoral and healthy regions.
7. Place a thin layer of OCT in a cryomold previously labeled and/or identified.
8. With sterile forceps, place 1 fragment into the cryomold with OCT. An adequate orientation in the cryomold is important to obtain optimal histological cuts in the cryostat.
9. Cover the lung tissue fragment with OCT.
10. Use forceps or a pipette tip to place the tissue properly oriented and remove any air bubbles.

11. Cool down isopentane holding its container over liquid nitrogen or dry ice. Isopentane is cool enough when it appears like pearls and the solution gets a white and dense appearance.
12. Place the cryomold into pre-cooled (-50°C) isopentane (in histobath or liquid nitrogen bath) with the aid of forceps. Avoid contact between isopentane and the OCT, to prevent air bubbles and lung tissue tearing.
13. Once frozen, transfer the cryomold to a -80°C freezer. Use liquid nitrogen or dry ice for transportation.
14. Register the stored sample location in the Pulmonary Biobank Consortium's software.

Protocols for tissue preservation are available in PNT 8.2.002, 8.2.003, 8.2.004, 8.2.005 y 8.2.006.

8. APPLICABLE REFERENCES, REGULATION AND GUIDELINES

1. Declaration of Helsinki. <http://ohsr.od.nih.gov/helsinki.php3>
<http://www.wma.net/e/policy/b3.htm>
2. Tri-Council Policy Statement; Ethical Conduct for Research Involving Humans; Medical Research Council of Canada; Natural Sciences and Engineering Council of Canada; Social Sciences and Humanities Research Council of Canada, August 1998.
<http://www.pre.ethics.gc.ca/english/policystatement/policystatement.cfm>
3. Human Tissue and Biological Samples for use in Research. Operational and Ethical Guidelines. Medical Research Council Ethics Series. http://www.mrc.ac.uk/pdf-tissue_guide_fin.pdf
4. Best Practices for Repositories I. Collection, Storage and Retrieval of Human Biological Materials for Research. International Society for Biological and Environmental Repositories (ISBER). <http://www.isber.org>
5. National Bioethics Advisory Commission: Research involving human biological materials: Ethical issues and policy guidance, Vol. I: Report and recommendations of the National Bioethics Advisory Committee. August 1999.
<http://bioethics.georgetown.edu/nbac/hbm.pdf>
6. US National Biospecimen Network Blueprint
http://www.ndoc.org/about_ndc/reports/NBN_comment.asp
7. Jewell, S. et al. 2002, Analysis of the Molecular Quality of Human Tissues, an experience from the Cooperative Human Tissue Network. Am. J. Clin.Pathol.118:733-741.
8. Guideline – Fresh Tissue Working Group of BIG and NCI breast cancer Cooperative Groups http://ctep.cancer.gov/forms/guidelines_fresh_tissue.pdf
9. SOP-PRO-Tissue-freezing.02, Oct 3, 2006. Freezing of Prostate Tissues. Procure
10. Quebec Prostate Cancer Biobank.